


How to install DinoCapture on Vista x64 OS


Step.1

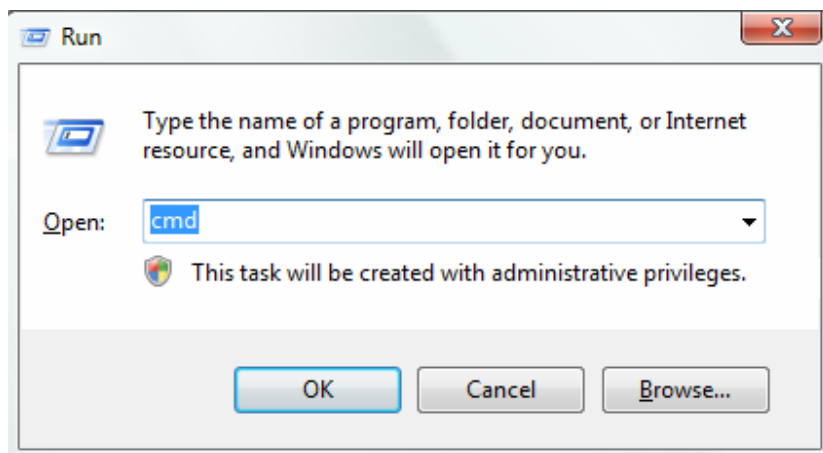



Step.2




Step.3

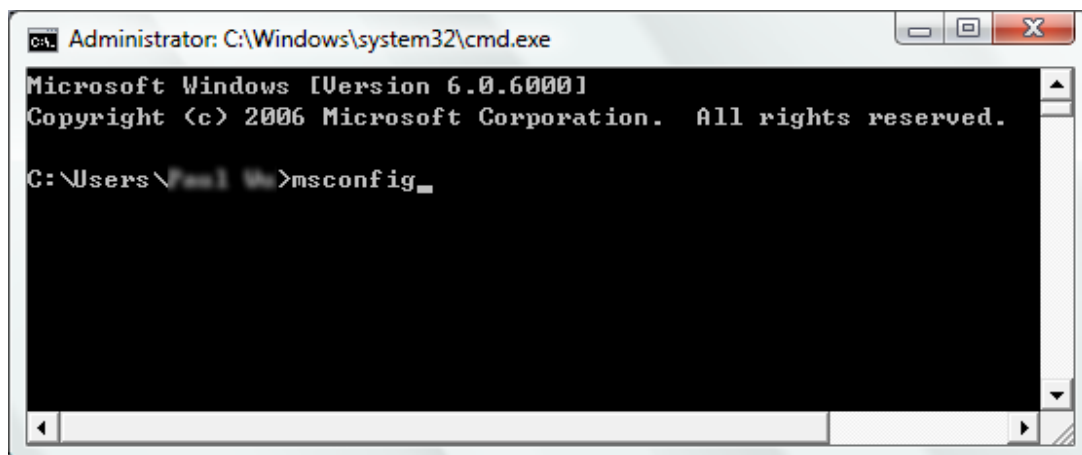


Step.4

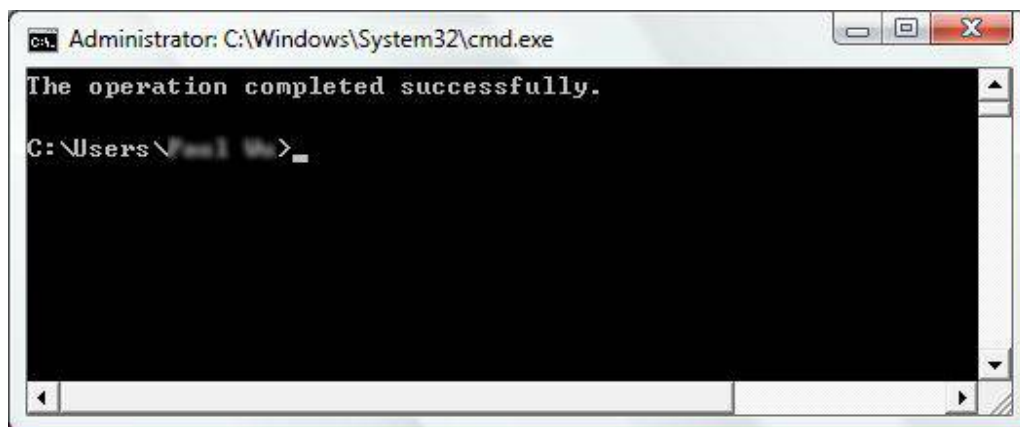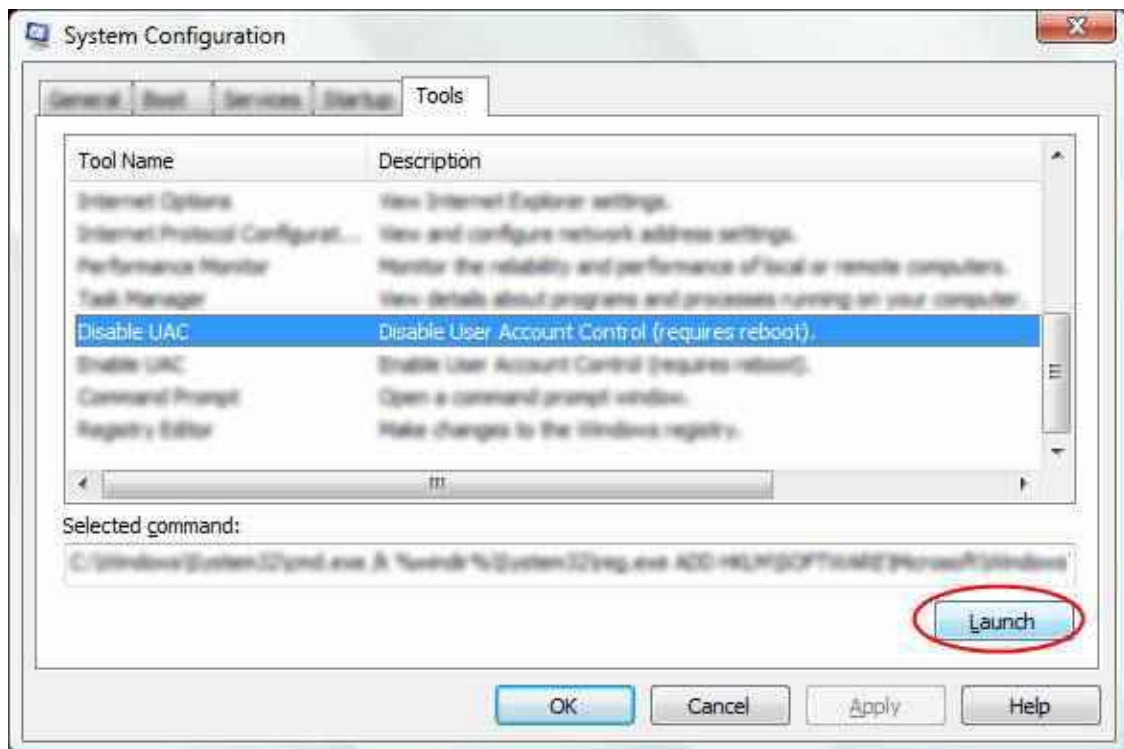

Step.5

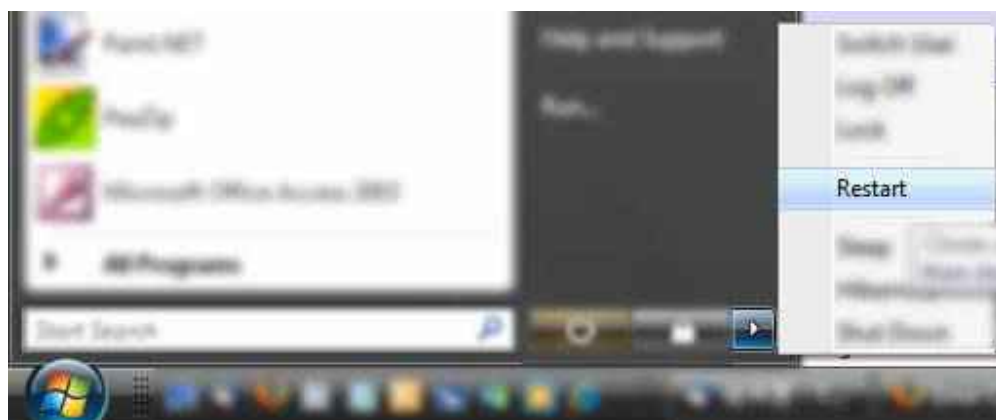

Step.6

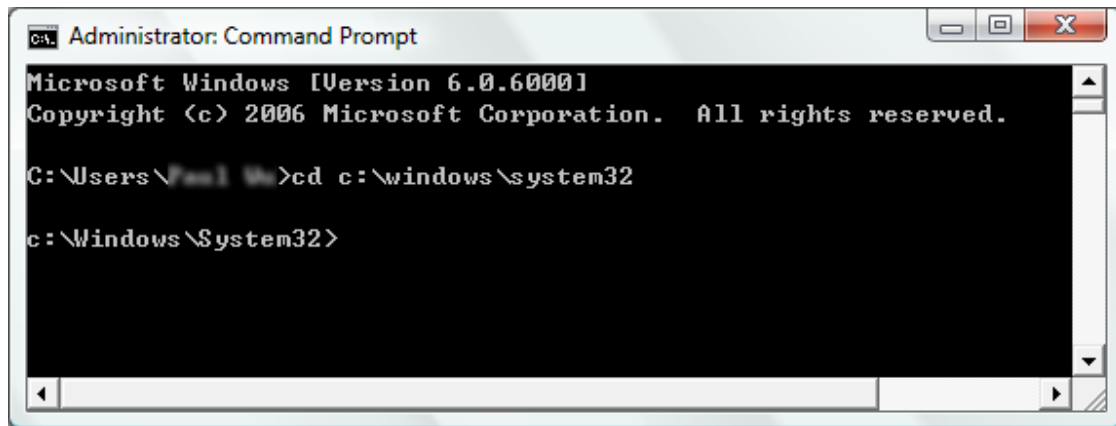


```
Administrator: Command Prompt
Microsoft Windows [Version 6.0.6000]
Copyright (c) 2006 Microsoft Corporation. All rights reserved.

C:\Users\Paul W>cd c:\windows\system32

c:\Windows\System32>
```

Step.7

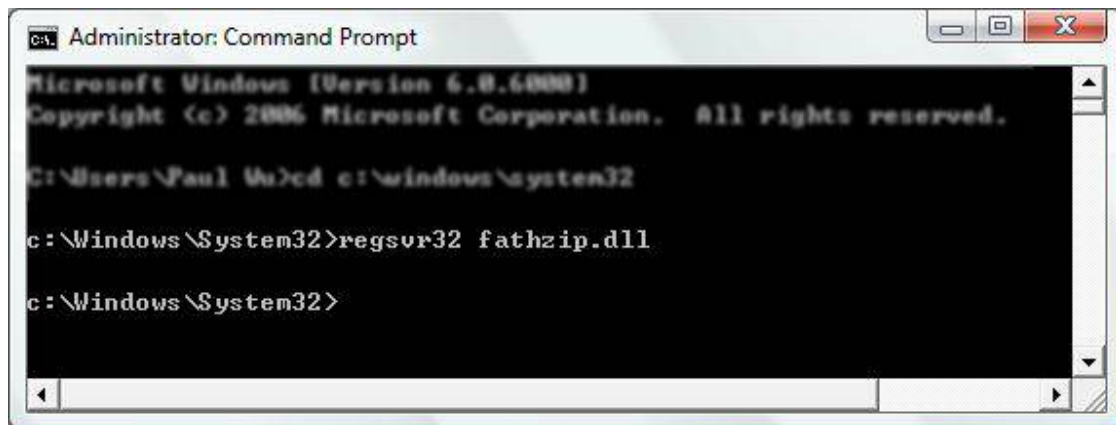


```
Administrator: Command Prompt
Microsoft Windows [Version 6.0.6000]
Copyright (c) 2006 Microsoft Corporation. All rights reserved.

C:\Users\Paul W>cd c:\windows\system32

c:\Windows\System32>regsvr32 fathzip.dll

c:\Windows\System32>
```

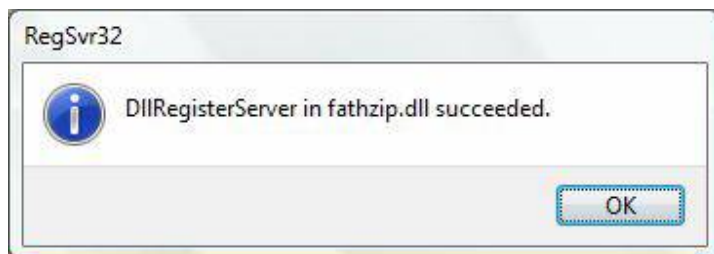

Step.8

Install the DinoCapture program.

(If the Dinocapture had installed before, please uninstall the DinoCapture first)